

How To Replace The Daylight Plate On The Refractometer

Refractometer RHCN-200ATC


Catalogue Number: 1841



Firstly, it is important to know which model of Refractometer you have purchased. At Vetlab Supplies Ltd we sell the **RHCN-200ACT** portable refractometer and can supply the correct replacement **daylight plate** for this model.

Step 1

Carefully remove the two screws using a small flat-headed screwdriver.

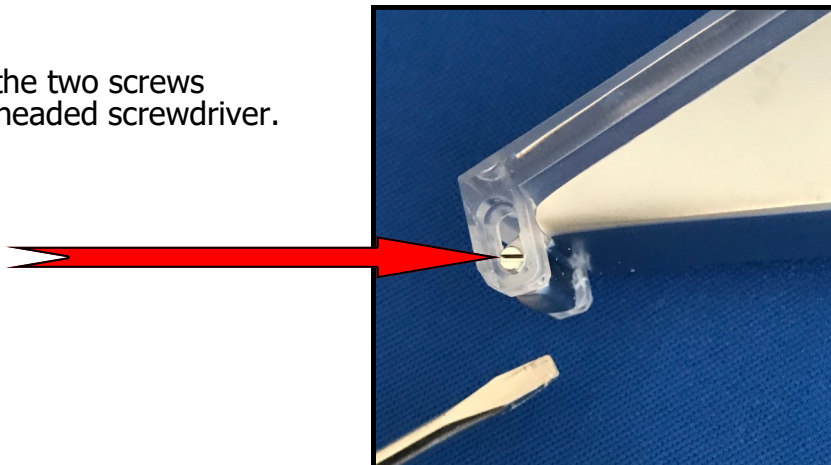

Step 2

Once you have removed the screws and the old daylight plate, carefully place the new daylight plate into the correct position and slowly tighten-up the screws.



The replacement daylight plate is supplied with two spare screws.

Refractometer RHCN-200ATC Daylight Plate

Catalogue Number: 1845

Please do not hesitate to contact us if you have any questions or require further assistance:

Telephone: 01798 874567

Email: info@vetlabsupplies.co.uk

We also supply: Veterinary Diagnostic Test Kits, Veterinary Equipment, Laboratory Consumables and Much More. Visit our website www.vetlabsupplies.co.uk