

MiniSpin NG Quick Guide to Changing Rotors

Please Follow Instructions Carefully - **NO TOOLS REQUIRED!**

USING/CHANGING THE 4-PLACE BLACK TUBE ROTOR



STEP 1-

Attaching Rotor

Hold the Rotor over the black clips. Line up correctly, push down and **CLICK** on securely.




STEP 2- Removing Rotor

Firstly, remove the Tube-Holders. Place the fingers of both hands under the rotor then using both thumbs squeeze in the **TOP** of the black clips and lift rotor off **VERTICALLY!**



IMPORTANT! DO NOT USE TOOLS OF ANY DESCRIPTION TO REMOVE ROTORS... SQUEEZE & LIFT!

USING/CHANGING THE MICROHAEMATOCRIT ROTOR

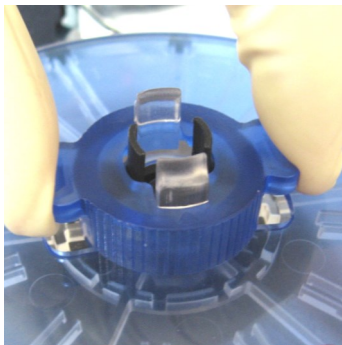

STEP 1 - Attaching Rotor


Firstly—REMOVE LID

Start by holding the larger wings with your finger and thumb.




STEP 1.1- Removing Lid

Using your other finger and thumb squeeze in/pinch in the other clips simultaneously and lift off the lid.

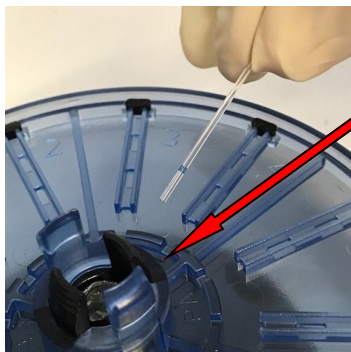


STEP 2 - Setting Up

Line up the **BASE** of the rotor with the black markers and push down until the rotor clicks on. Load your samples – **Sealant end outwards.**

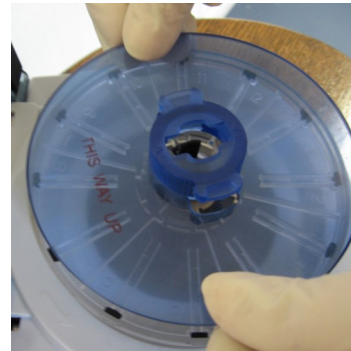

STEP 3 - Replacing Lid

Always replace the rotor lid for PCV's! Line it up with the black markers and **CLICK** down. Set your Timer and switch on.

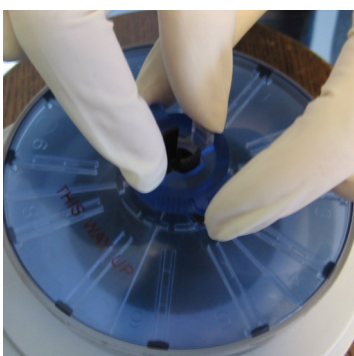

STEP 4 - Removing Lid

After spinning down, remove the rotor lid. **See Step 1**, leaving the capillaries in the base section. Remove the capillaries and read.



STEP 5 - Removing Base

To remove the base place the fingers of both hands under the rotor using your thumbs to squeeze/pinch the **TOP** of the black clips inwards and lift the rotor off vertically.

SPINNING TIMES: PLEASE ALWAYS USE A TIMER!

PLASMA/SERUM 2 Minutes Minimum: MICROHAEMATOCRIT 3 Minutes Minimum: URINE 20 Seconds:

Please contact technical services on: 01798 874871 if you need any assistance. (16.08.17)