

VetSpin® Duo Quick User Guide

Please follow instructions carefully



START UP

A. Prepare VetSpin

- Switch on at back of unit
 - Press **STOP/LID** key to OPEN LID
- DO NOT TOUCH THE TOP OF THE LID WITH FINGERS/HANDS WHILE PRESSING THE **STOP/LID** KEY



B. Remove Rotor Cover

(if the Rotor Cover is fitted) Unscrew clear plastic rotor cover and remove ready to load samples. Ensure all **TUBE HOLDERS** are in place. **Never run with missing TUBE HOLDERS!**

The Rotor Nut secures the rotor onto the spindle and must be tightly in place at all times! **Use Your Allen Key To Tighten.**

Serum/Plasma
Load Tubes and Remember to **BALANCE!**
MAX 2ml Tubes




TUBES
Never remove tube holders to try and accommodate incorrect sized tubes!




PCV/(MHct) 75mm
Load Capillaries – sealant end outwards & fit the **ROTOR COVER!**



Tubes & PCV's
Load Tubes and Remember to **BALANCE & fit the ROTOR COVER!**

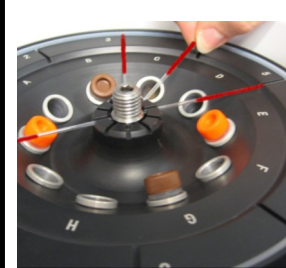

Urine
Load Tubes and Remember to **BALANCE!**

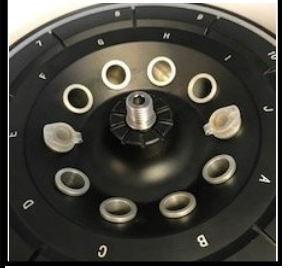

1. Press **YELLOW** Program Button
2. Select Program Required

Serum/Plasma Blood Tubes
Pre-programmed 2 Mins FAST SPEED

MAX SIZE 2ml PLASTIC only - Never remove tube holders to accommodate incorrect sized tubes, this could cause catastrophic damage to the centrifuge. No Glass - No Vacutainers.



PCV (MHct) - 75mm length
Pre-programmed 5 Mins FAST SPEED OR - 2 Mins Can Be Enough

Use also for Combination Spinning
(Blood Tubes & PCV/MHct Together)

Urine Sediment - Pre-programmed 2 Mins SLOW SPEED



- Once you have loaded your tubes correctly, close the lid and press the **GREEN START KEY.**
- Stay with the centrifuge until it has reached full speed in case it needs to be turned off! **RED STOP KEY!**
- **TURN OFF IMMEDIATELY if there is any wobble, shaking or unusual noise which will be caused by incorrect balancing. Incorrect balancing can seriously damage the centrifuge.**
- When the run has finished the centrifuge will BEEP, **3 Triple Beeps.** Press **RED** button to open lid.
- Remove the samples and leave the lid open so it can air, cool down and remain dry.

This is a Quick User Guide Only - Please use the full User Manual for detailed information.
Contact: Technical Support on: 01798 874567 if you have any queries or need assistance. VetSpin Dec 2020