

# VetSpin® Duo Quick User Guide

Please Follow Instructions Carefully



## START UP

### A. Prepare VetSpin

- Switch on at back of unit
- Press **STOP/LID** key to OPEN LID

DO NOT TOUCH THE TOP OF THE LID WITH FINGERS/HANDS WHILE PRESSING THE **STOP/LID** KEY



### B. Remove Rotor Cover

(if the Rotor Cover is fitted) Unscrew clear plastic rotor cover and remove ready to load samples. Ensure all **TUBE HOLDERS** are in place. Never run with missing **TUBE HOLDERS!**

**The Rotor Nut** secures the rotor onto the spindle and must be tightly in place at all times! **Use Your Allen Key To Tighten.**

### Serum/Plasma

Load Tubes and Remember to **BALANCE!**



### PCV/(MHct) 75mm

Load Capillaries – sealant end outwards & fit the **ROTOR COVER!**



### Tubes & PVC's

Load Tubes and Remember to **BALANCE & fit the ROTOR COVER!**



### Urine

Load Tubes and Remember to **BALANCE!**



1. Select Program Required
2. Press **YELLOW** Program Button

### Serum/Plasma

*Pre-programmed 2 Mins FAST SPEED*

### PCV (MHct) - 75mm length

*Pre-programmed 5 Mins FAST SPEED*

**Use also for Combination Spinning**  
(Blood Tubes & PCV/MHct Together)

### Urine Sediment

*Pre-programmed 2 Mins SLOW SPEED*

- Once you have loaded your tubes correctly, close the lid and press the **GREEN START KEY.**
- Stay with the centrifuge until it has reached full speed in case it needs to be turned off! **RED STOP KEY!**
- **TURN OFF IMMEDIATELY if there is any wobble, shaking or unusual noise which will be caused by incorrect balancing. Incorrect balancing can seriously damage the centrifuge.**
- When the run has finished the centrifuge will BEEP, **3 Triple Beeps.** Press **RED** button to open lid.
- Remove the samples and leave the lid open so it can air and remain dry.

**This is a Quick User Guide Only - Please use the full User Manual for detailed information.**  
**Contact: Technical Support on: 01798 874871 if you have any queries or need assistance. New VetSpin 2019**