

# VetSpin® Micro Quick Guide to Changing Rotors

Please Follow Instructions Carefully - **NO TOOLS REQUIRED!**

## USING/CHANGING THE 4-PLACE BLACK TUBE ROTOR



### STEP 1-

#### Attaching Rotor

Hold the Rotor over the black clips. Line up correctly, push down and **CLICK** on securely.




### STEP 2- Removing Rotor

Firstly, remove the Tube-Holders. Place the fingers of both hands under the rotor then using both thumbs squeeze in the **TOP** of the black clips and lift rotor off **VERTICALLY!**



**IMPORTANT! DO NOT USE TOOLS OF ANY DESCRIPTION TO REMOVE ROTORS... SQUEEZE & LIFT!**

## USING/CHANGING THE MICROHAEMATOCRIT ROTOR

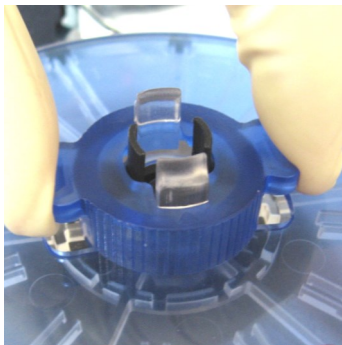

### STEP 1 - Attaching Rotor


#### **Firstly—REMOVE LID**

Start by holding the larger wings with your finger and thumb.




### STEP 1.1- Removing Lid

Using your other finger and thumb squeeze in/ pinch in the other clips simultaneously and lift off the lid.

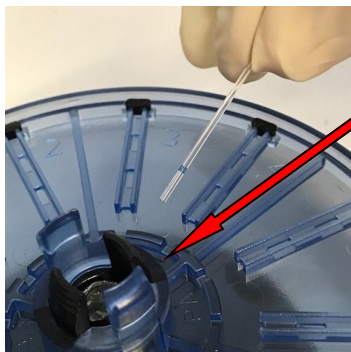


### STEP 2 - Setting Up

Line up the **BASE** of the rotor with the black markers and push down until the rotor clicks on. Load your samples – **Sealant end outwards.**

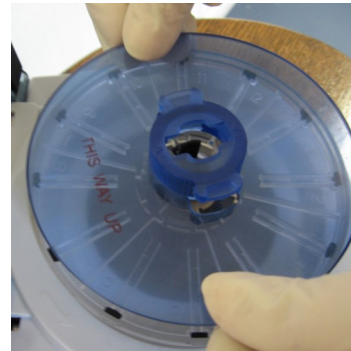

### STEP 3 - Replacing Lid

**Always** replace the rotor lid for PCV's! Line it up with the black markers and **CLICK** down. Set your Timer and switch on.

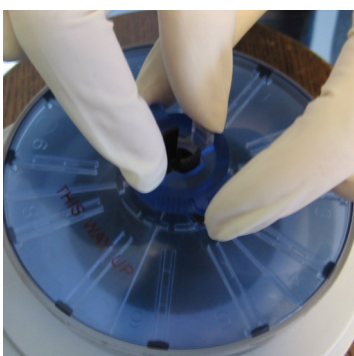

### STEP 4 - Removing Lid

After spinning down, remove the rotor lid. **See Step 1**, leaving the capillaries in the base section. Remove the capillaries and read.



### STEP 5 - Removing Base

To remove the base place the fingers of both hands under the rotor using your thumbs to squeeze/ pinch the **TOP** of the black clips inwards and lift the rotor off vertically.

**SPINNING TIMES: PLEASE ALWAYS USE A TIMER!**

**PLASMA/SERUM 2 Minutes Minimum: MICROHAEMATOCRIT 3 Minutes Minimum: URINE 20 Seconds:**

Please contact technical services on: 01798 874871 if you need any assistance. (March 2019)